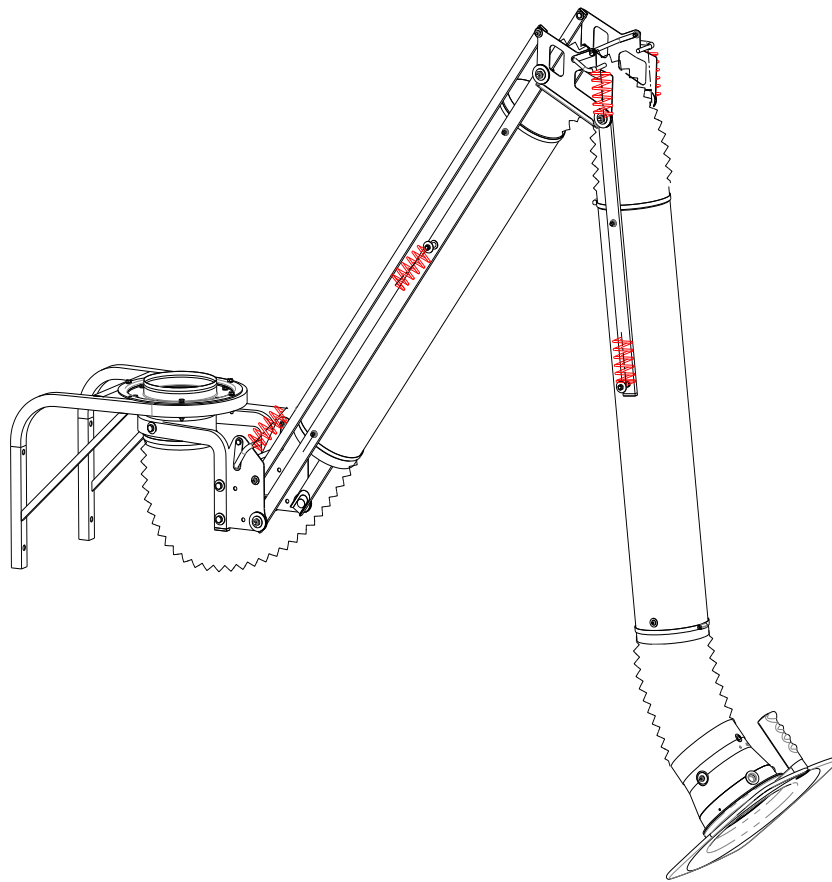
HRM160 series Suction Arms
Bras Aspirants série HRM160
Bracci Aspiranti serie HRM160
Absaugarme Serie HRM160

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EN

Reference standards:

- UNI EN 12100-2010 - Requirements for the design in machine safety
- UNI EN 13857:2020 - Safety distances to prevent danger zones being reached by the upper limbs

WARNINGS and METHODS OF USE	3
PACKAGING	5
INSTALLATION	7
Cabinet version suction arm assembly.....	7
Standing version suction arm assembly.....	11
MAINTENANCE	12

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WARNINGS and METHODS OF USE

- ! Please read carefully the regulations described below before starting to assemble the arm.
- ! **Before starting to unpack, install or perform maintenance, equip yourself with suitable personal protective equipment, i.e. gloves, shoes and a safety helmet, and the following tools: cutter, 2 x 10 mm Allen wrenches and 2 x 13 mm Allen wrenches.**
- ! The arm must only be handled by using the specific handle on the hood.
- ! The position of the hood must be coordinated with the suction flow rate and in any case no more than 30 cm from the source of pollution.
- ! Pay utmost attention when handling the arm so as not to accidentally bump into objects and/or people.
- ! Under no circumstances must the liquid and/or gas over 90°C (except for special models for high temperatures), explosive compounds (e.g. titanium or magnesium powder, etc.), hot and/or inflamed objects be drawn or even draw near open flames.
- ! Under no circumstances is suction to be performed with a vacuum exceeding 1600 Pa (except for special high-vacuum models).

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(1) ARM BODY WITH ROTATING BASE (EN)
 (2) BRACKET FOR WALL MOUNTING
 (3) FITTING
 (4) HOOD
 (5) BAG WITH INSTALLATION KIT AND MANUAL

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(*)=Standing Version obtained by transforming the cabinet version.

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EN

PACKAGING

Fig. A shows the components inside the packaging with the weights and dimensions shown in Tab. 1.

An adhesive label indicating the following information is found on the outer part of the packaging:

- Sender, recipient, delivery address, type of goods contained and other information required by the carrier.
-  Before handling the packed product from the delivery vehicle, check that the packaging is entirely intact so as to prevent any components from leaking and/or falling out.
-  When unpacking, move the removed material away from the operating areas so as to eliminate the risk of the operator stumbling over it.
-  Under no circumstances must the end components of the package (cardboard, polyethylene, metal) be disposed of in the environment.
-  It is prescribed for disposal to be implemented by using the regular waste disposal channels according to local regulations.

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EN

ASSEMBLY ON WALL MOUNTED BRACKET

- Fig. B – Secure the bracket to the wall according to the values specified in the figure.
Before mounting the bracket, make sure that the wall is sufficiently flat and solid for the arm.
 While installing the bracket, verify that it does not deform and that it is horizontal.
An imprecise assembly can cause rotary difficulties or in an extreme case, the rotation to block.
- Fig. C – ALT. 1 and 2: Assemble the fitting or the fan by removing the 4 M6x40 screws and then reassemble them (in the case of a fan, the nuts are disposable).
- Raise the arm until the flange (1) of the bracket base is rested, while inserting the seal (2); then secure it using the 6 M6x16 screws supplied in the bag.

 Risk of shearing the fingers if the base rotates inadvertently.

ASSEMBLY ON JIB EXTENSION

- Fig. C (ALT.3) - Proceed by installing the arm on the jib extension as described in point 1.2 above. Verify that the configuration of the holes on the jib extension is compatible with those of the base (view A-A).

 Risk of shearing the fingers if the base rotates inadvertently.

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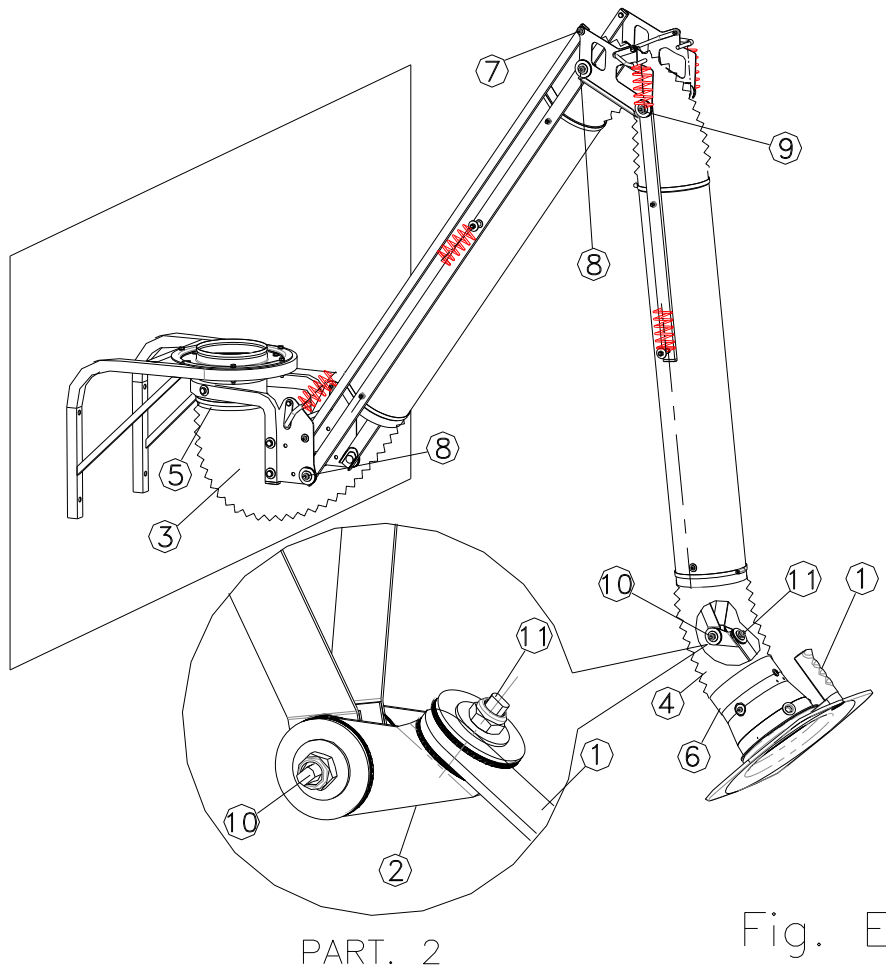
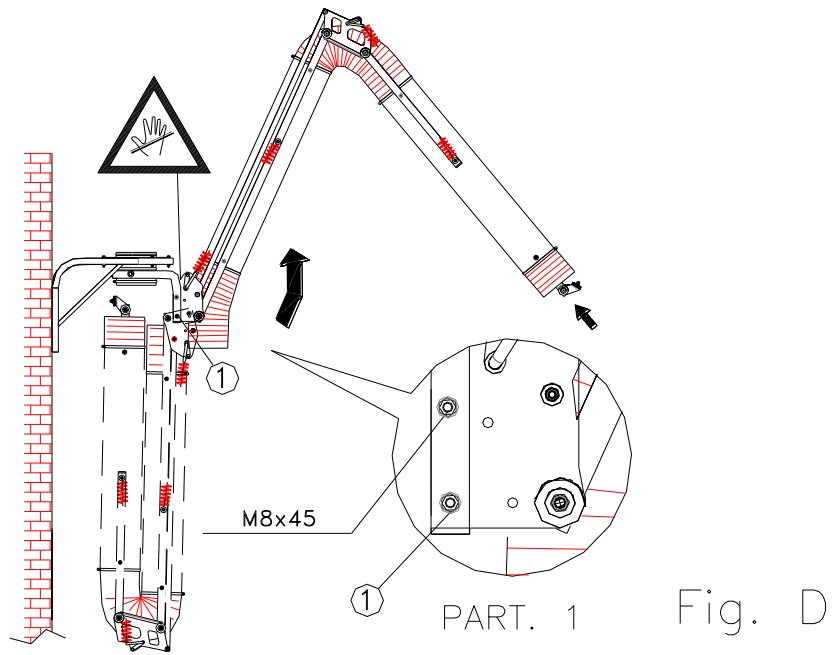
INSTALLATION - Assembly in cabinet version

EN

- 1) Fig. D - Raise the arm body by rotating it upwards until the 2 M8x45 screws can be inserted (DET. 1).
 Risk of shearing fingers between the base and the plates of the arm.
- 2) Fig. E - Assemble the hood (1) on the double bracket (2) according to the assembly diagram shown in the figure (DET. 2). Adjust the arm by tightening or loosening the joint nuts (7), (8), (9), (10) and (11) by 1/8 turn. Finally secure the flexible hoses (3-4) to the base and to the hood using the clamp (5) and the rubber collar (6).

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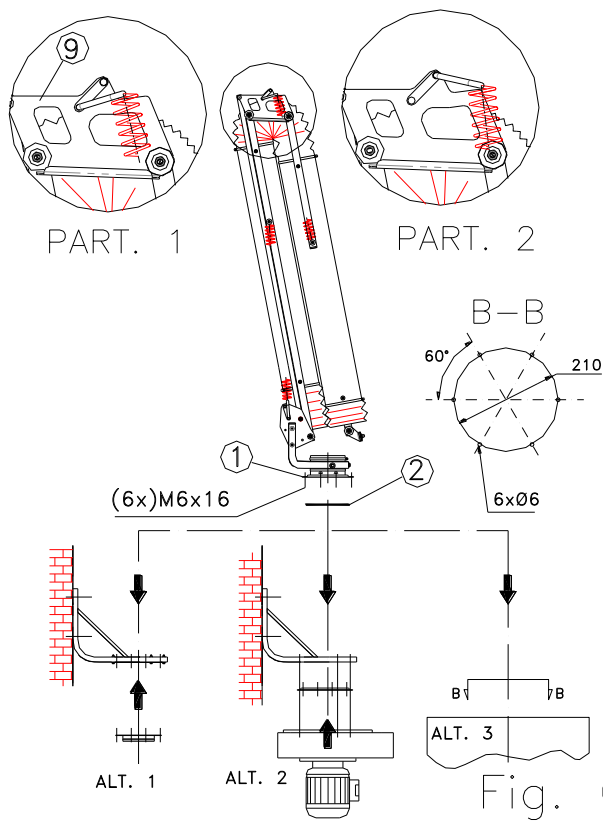
Fig. G

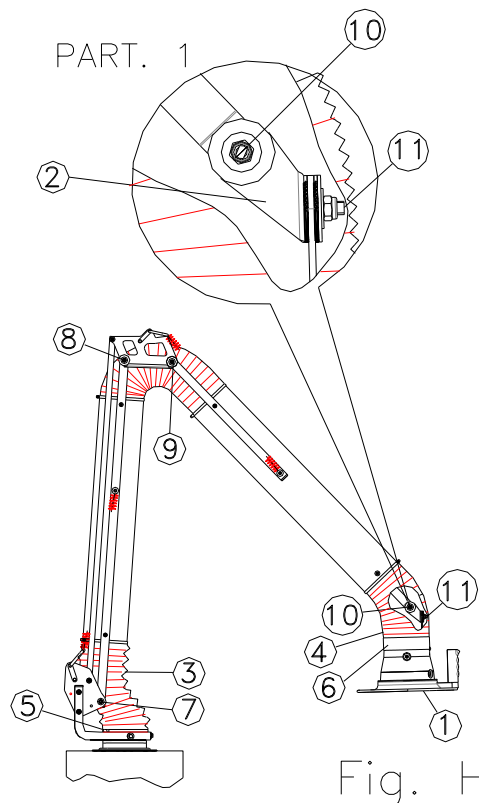
Fig. H

INSTALLATION - Transformation from the Cabinet to Standing version (Only for 2 and 3m; request for 4m.)

EN

- 1) Fig. F - Remove the base (1) from the arm by unscrewing the 2 M8x45 screws to reassemble it in the new position (Det.1-2)
- ⚠ CAUTION! Risk of shearing fingers between the tubes of the base and the plates of the arm body.
- 2) Fig. F - Loosen the nut (3) and release the spring (4) from the pin. Slide the spring coupling hook (6) downwards towards the lower part of the slot (7) (Det. 3) after having loosened the screw (5) which blocks the cover-slot plate. Insert the spacer (8) by coupling the spring (Det. 4). Repeat for each spring.
- 3) Fig. G - Repeat the sliding steps of the hook also on the trapeze (9) but without inserting the spacer (Det.1-2).

ASSEMBLY ON TROLLEY

Fig. G - Prepare the holes on the trolley as indicated in Sec. B-B. Assemble the arm by placing the seal (2) and using the 6 M6x16 screws supplied in the installation kit.

Make sure that the trolley structure is sufficiently robust and stable to prevent it from tipping over.

ASSEMBLY ON BRACKET Fig. H - Refer to the instructions for the cabinet version on page 9.

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MAINTENANCE

Maintenance must be performed by qualified and skilled personnel by reading this manual.

EN

POS.	PART	DUE DATE	MAINTENANCE
1)	FRICTION	If the arm is unstable	<ul style="list-style-type: none"> Adjust the friction of the joints by loosening or tightening the blocking nuts (Fig. G, Fig. L).
2)	BASE SEAL	12 months	<ul style="list-style-type: none"> Clean any debris from the seals and lubricate everything before assembly (Fig. D, Fig. E, Fig. J, Fig. K).
3)	FLEXIBLE HOSES	12 months	<ul style="list-style-type: none"> Check that the flexible hose is intact and the clamps and/or rubber collars are sealed; if necessary, replace them.

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