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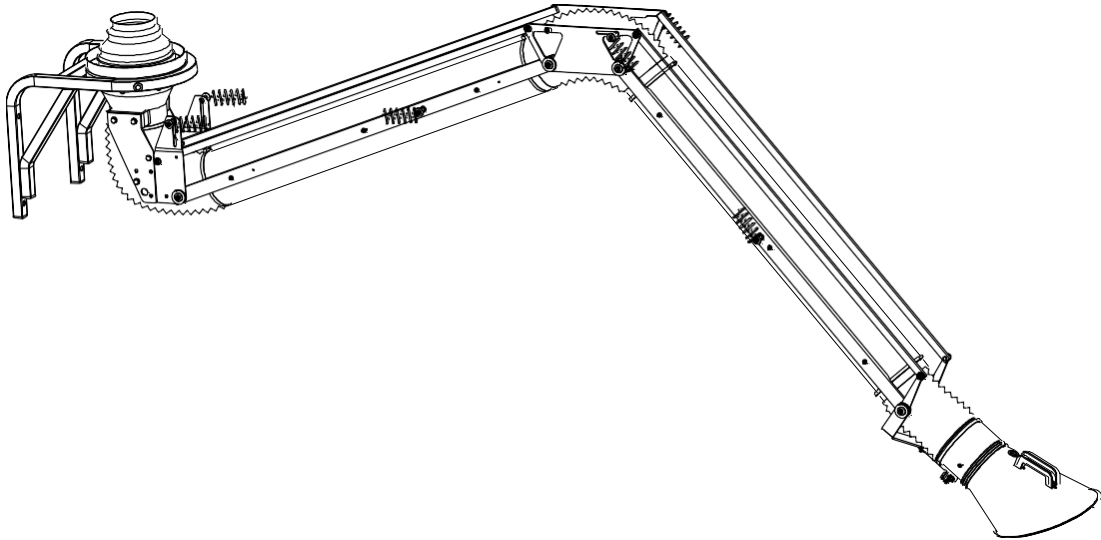
EN

Reference standards:

- UNI EN 12100-2010 - Requirements for the design in machine safety
- UNI EN 13857:2008 - Safety distances to prevent danger zones being reached by the upper limbs.

1	INDEX	2
2	WARNINGS and METHODS OF USE	3
3	INSTALLATION.....	4
3.1	Hanging version suction arm assembly	4
3.2	Standing version suction arm assembly.....	8
4	MAINTENANCE.....	12
5	PACKAGING.....	13

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WARNINGS and METHODS OF USE

EN

- ! Please read carefully the regulations described below before starting to assemble the arm.
- ! Before starting to unpack, install or perform maintenance, equip yourself with suitable personal protective equipment, i.e. gloves, shoes and a safety helmet, and the following tools: cutter, 2 x 10 mm and 2 x 13 mm Allen wrenches.
- ! The arm must only be handled by using the specific handle on the hood.
- ! The position of the hood must be coordinated with the suction flow rate and in any case not further than 30 cm from the source of pollution.
- ⚠ Pay utmost attention when handling the arm so as not to accidentally bump into objects and/or people.
- Under no circumstances must the liquid and/or gas over 90°C (except for special models for high temperatures), explosive compounds (e.g. titanium or magnesium powder, etc.), hot and/or inflamed objects be drawn or even draw near open flames.
- Avoid to apply to the suction arm a vacuum exceeding 1600Pa (except for special high-vacuum models).

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INSTALLATION
 INSTALLATION
 INSTALLAZIONE
 INSTALLATION

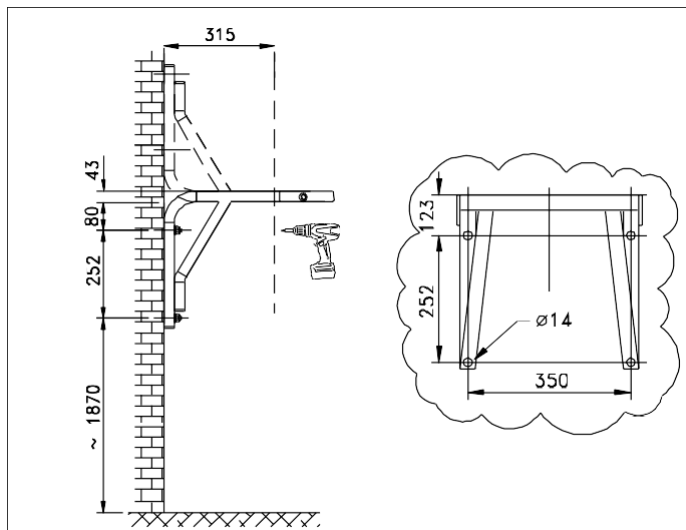
Fig./Abb. A

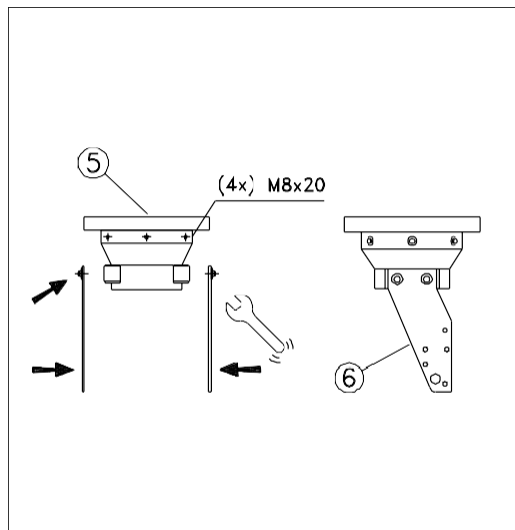
Fig./ Abb. A1

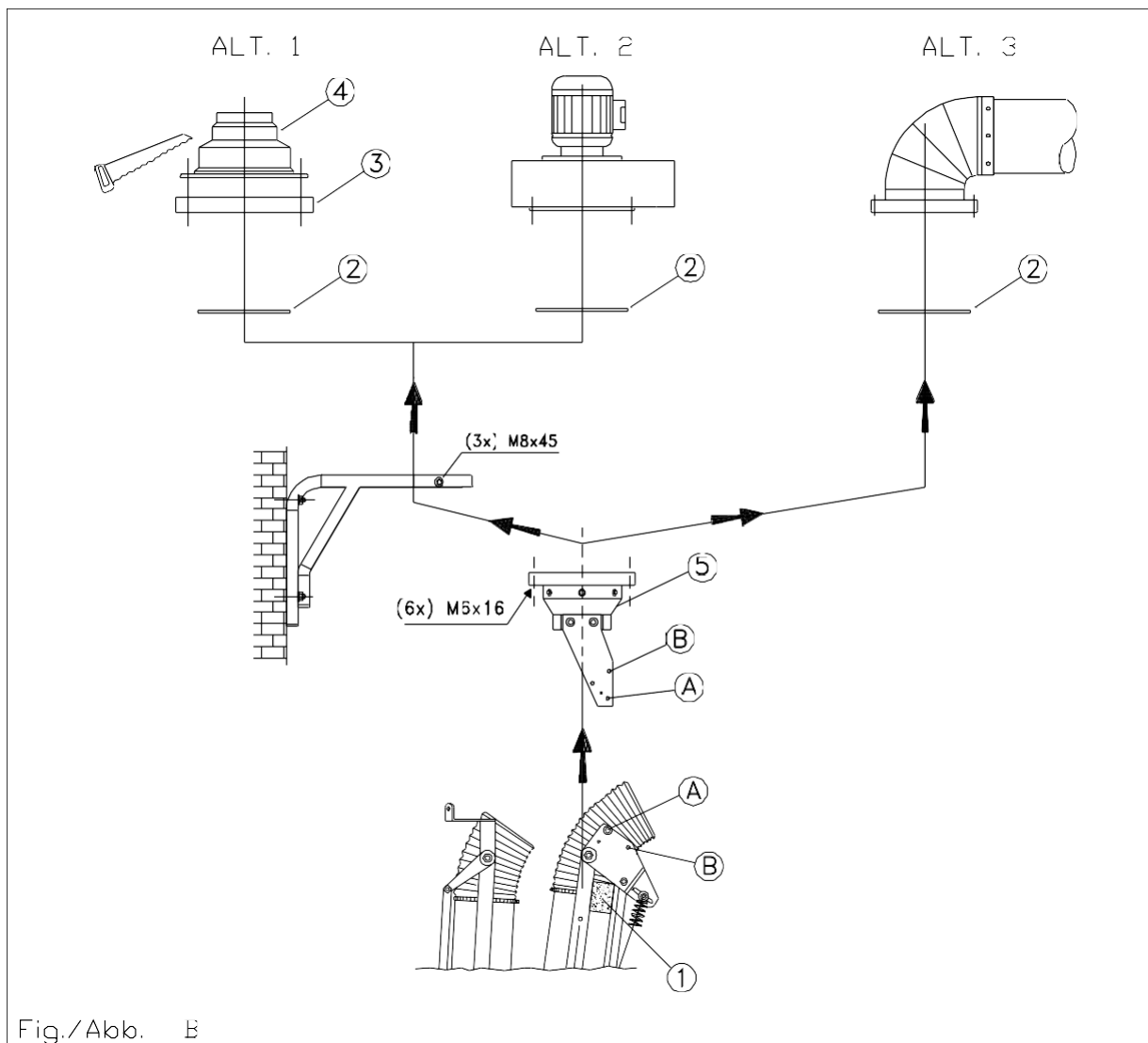
Fig./Abb. B

**INSTALLATION
INSTALLATION
INSTALLAZIONE
INSTALLATION**

ASSEMBLY OF THE HANGING MODEL

EN

- ! Fig. A - Before the installation of the wall-bracket, make sure that the wall is sufficiently flat and sturdy to provide proper bearing. During the installation make sure that the bracket does not become deformed and that it is positioned in perfectly horizontal alignment. Inaccurate assembly can cause rotation problems or the block of the rotation itself. Fix the wall-bracket to the height shown in the figure.
Fig. A1 – Assemble the 2 mounting brackets (6) on the revolving base plate (5) with the 4 M8x20 screws
- ! Fig. B (ALT. 1 and 2) - Insert the revolving base plate (5) in the wall bracket and fix it in place with the 3 M8x45 screws included in the installation kit bag. Assemble the counter flange (3) and the multi-diameter (4) cut at the appropriate diameter (ALT. 1) or the fan (ALT.2) using the 6 M6x12 screws included. Do not forget to insert the gasket (2).
- ! Fig. B (ALT. 3) – JIB EXTENSION MOUNTED VERSION - Assemble the revolving base plate on the jib extension previously fixed using the 6 M6x12 screws included in the installation kit bag. Do not forget to insert the gasket (2).
- Extract the arm body from the package and draw it near to the previously installed base. Do not remove the spacers (1) until the assembly is complete.

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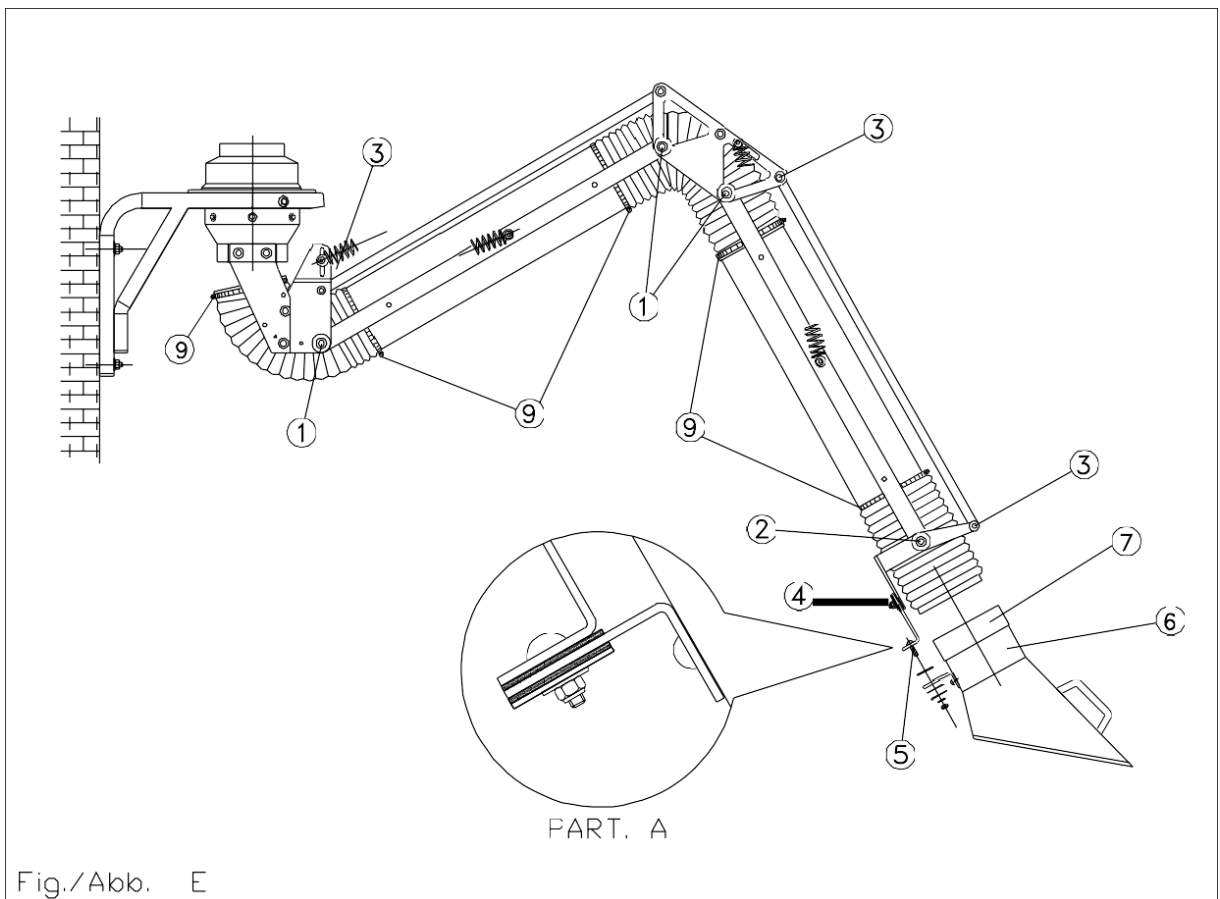
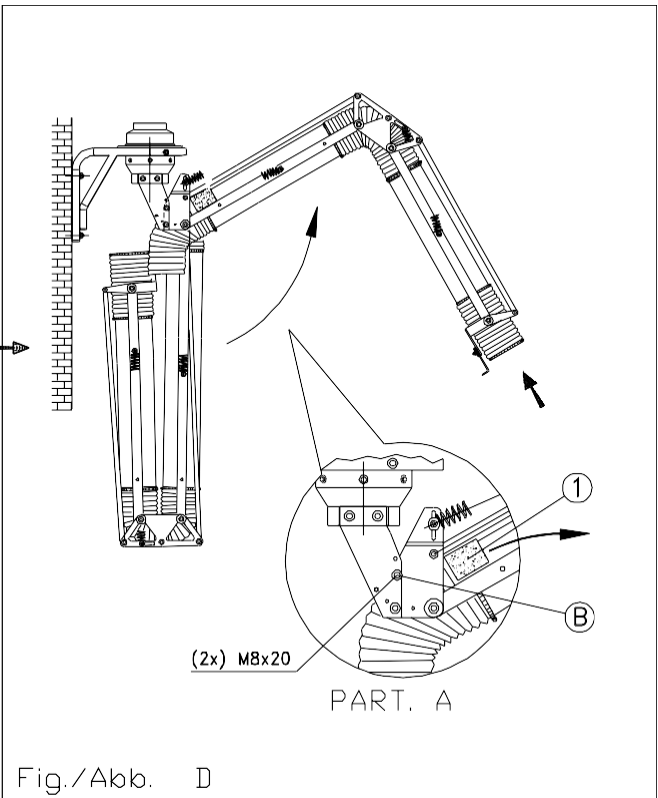
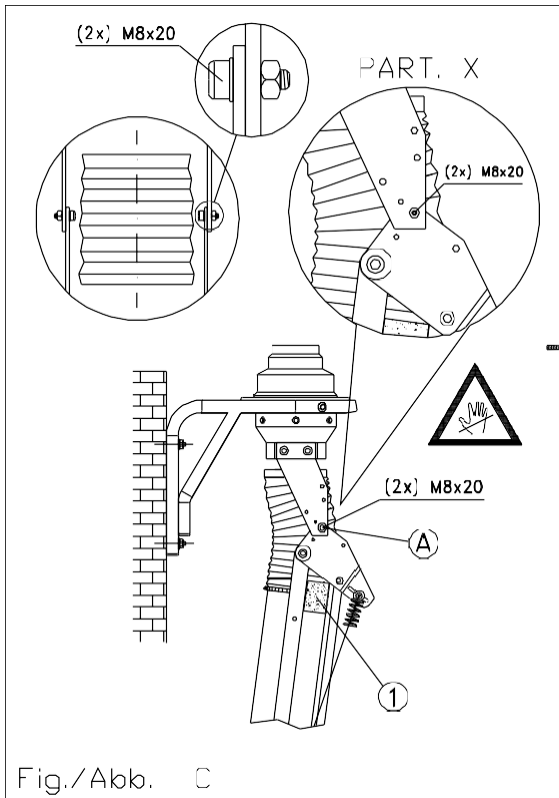
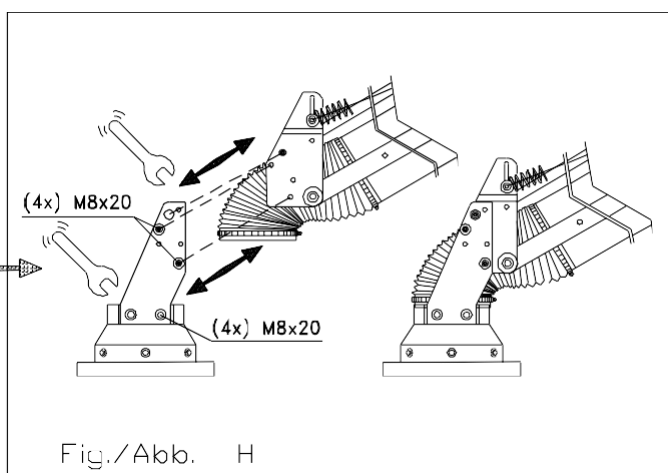
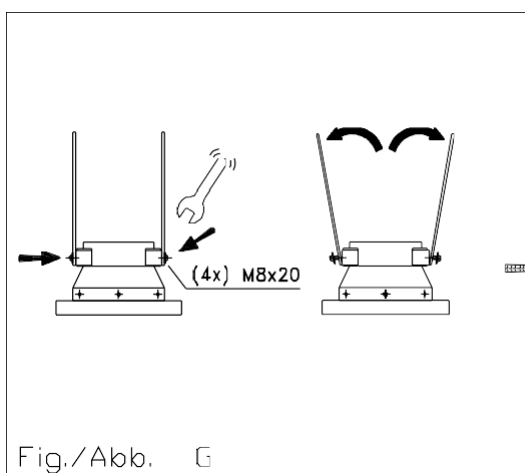
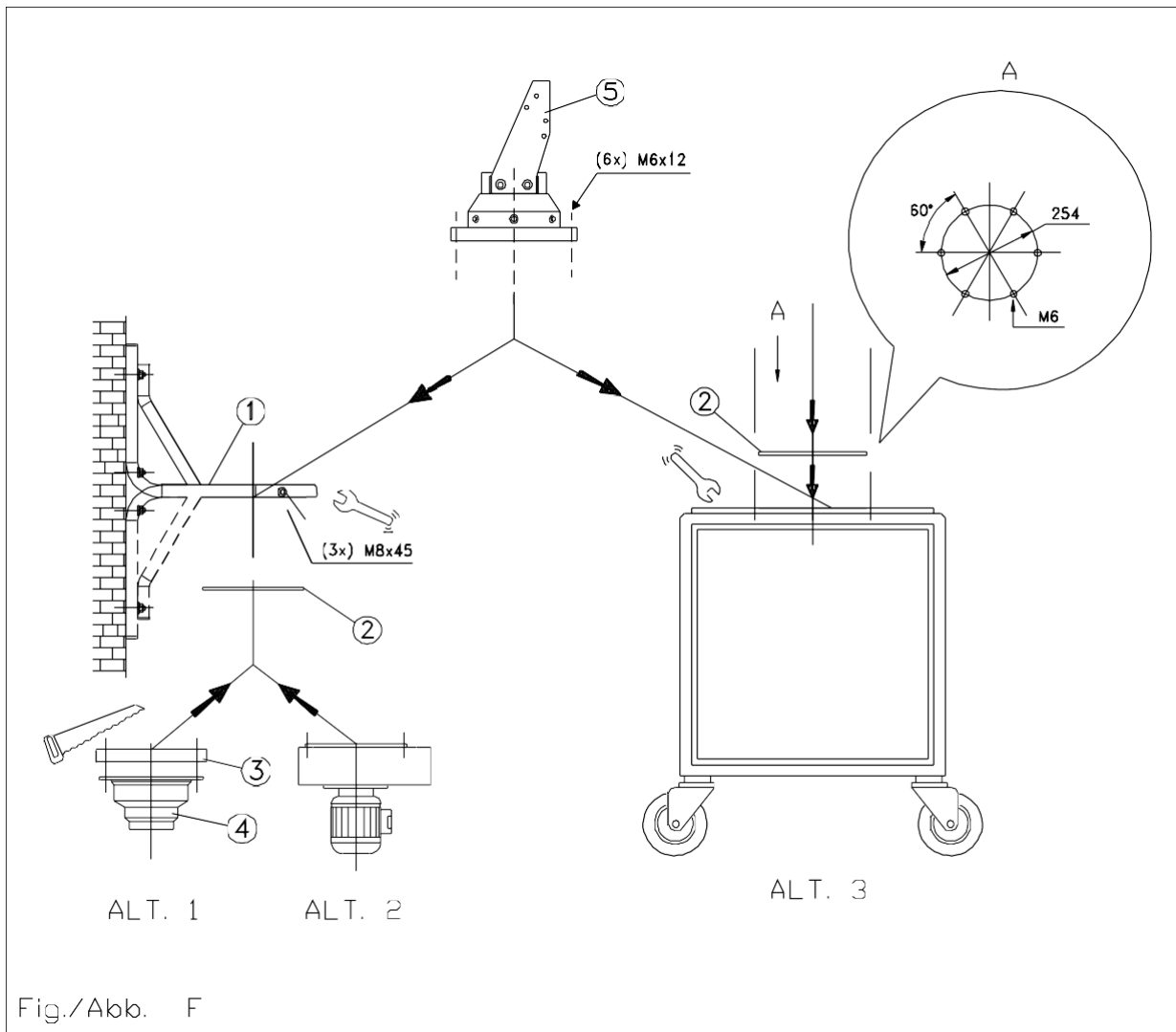
INSTALLATION
 INSTALLATION
 INSTALLAZIONE
 INSTALLATION

**INSTALLATION
INSTALLATION
INSTALLAZIONE
INSTALLATION**

STANDING VERSION

EN

- ! Fig. F (ALT. 1 e 2) - Fix the wall-bracket to the wall and assemble the revolving base plate as required on page 5 (Fig. A e A1). Insert the revolving base plate in the wall bracket (Fig. F) and fix it in place with the 3 M8x45 screws included in the installation kit bag. Assemble the counter flange and the multi-diameter cut at the appropriate diameter (ALT. 1) or the fan (ALT.2) using the 6 M6x12 screws included. Do not forget to insert the gasket (2).
Fig. F (ALT. 3) - Make sure that the carriage structure is sturdy and stable enough for the arm, then set up the drilling for the revolving base plate (Det. A) and fix it using the 6 M6x12 screws included. Do not forget to insert the gasket (2).
- ! Figs. G-H- Loosen the 4 M8x20 bolts of the base and extend the plates as shown (Fig. G); draw the arm body close, fixing it to the base with the 4 M8x20 bolts as shown in Fig. H and tighten the 4 base bolts again.
Fig. E (page 6) - Fix the hood (6) using the body square (5), the rubber collar (7) and clamps (9) as shown for the hanging version on page 7.

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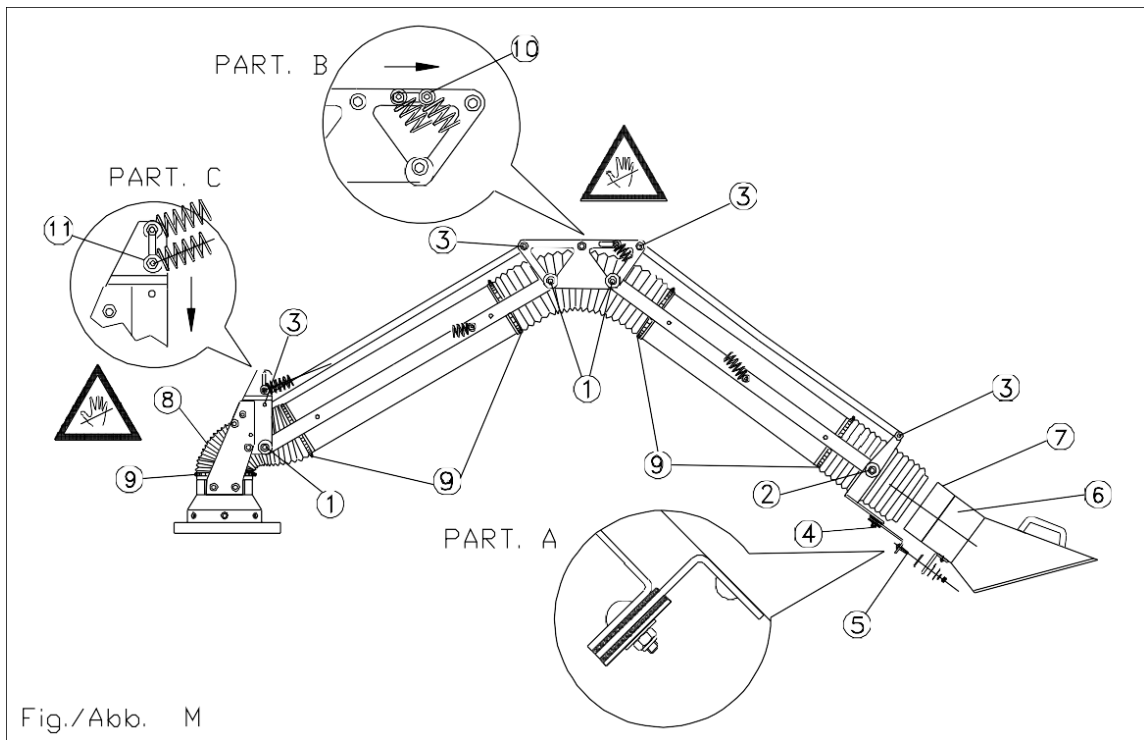
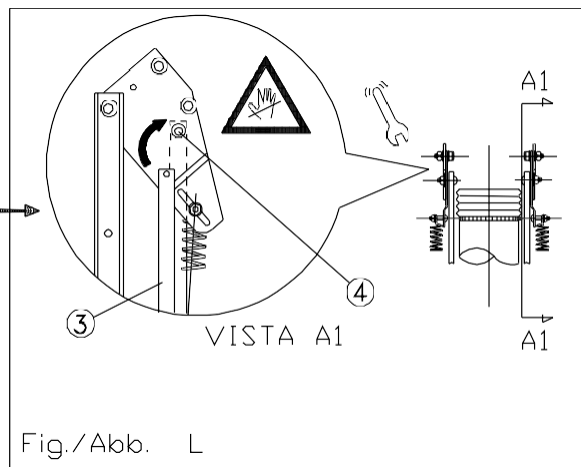
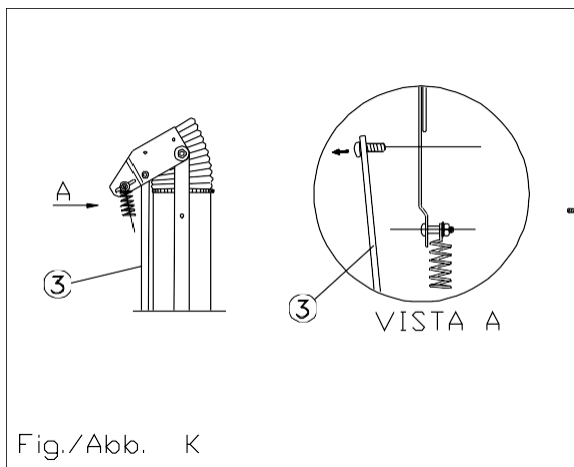
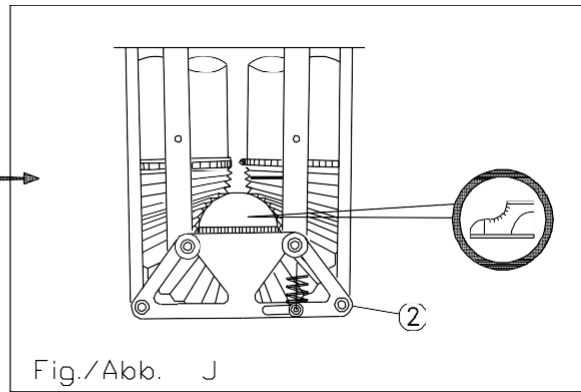
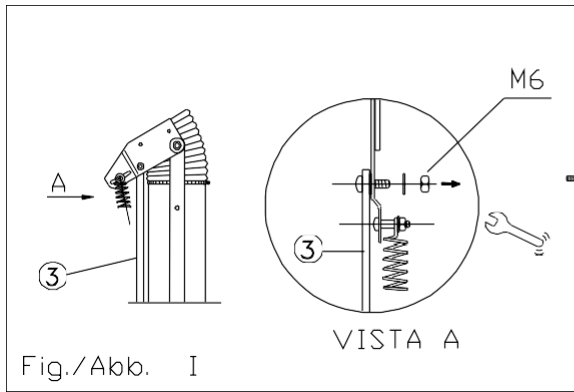
INSTALLATION
 INSTALLATION
 INSTALLAZIONE
 INSTALLATION

**INSTALLATION
INSTALLATION
INSTALLAZIONE
INSTALLATION**

CONVERTING HANGING VERSION INTO STANDING VERSION

EN

 Fig. I-J-K - Remove the M6 nut from the tie rod (3) (Fig. I). Place your foot on the piece (2) (Fig. J). Then, lift until you are able to remove the tie rod with its screw (Fig. K). Perform the same operations on the other side.

 Fig. L - Keeping your fingers in a safe position, let the piece (3) go gently until you can insert the screw in its new upper hole (4). Replace the M6 nut in the same order, then fasten. Perform the same operations on the other side.

Fig. F-H – At this point continue as per Fig. F to Fig. H on pages 8 and 9.

Fig. M - Loosen the nuts (10) by not more than one turn to advance both the anchoring of the springs on the trapezoidal arm piece (Part. B) (be careful not to loosen the nuts for more than one turn as the springs could be released) then re-tighten them. Loosen the fixing screws (11) of the springs by one turn (Part. C); to allow the fixing screws to slide down move the suction arm in its lowest position, then tighten the screws again. Cut the hose (8) which has to be connected with the base plate down by 30 cm. (~ 1 ft). Then, proceed as shown in the Fig. (E) on pages 6 and 7.

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**PACKAGING
EMBALLAGE
IMBALLAGGIO
VERPACKUNG**

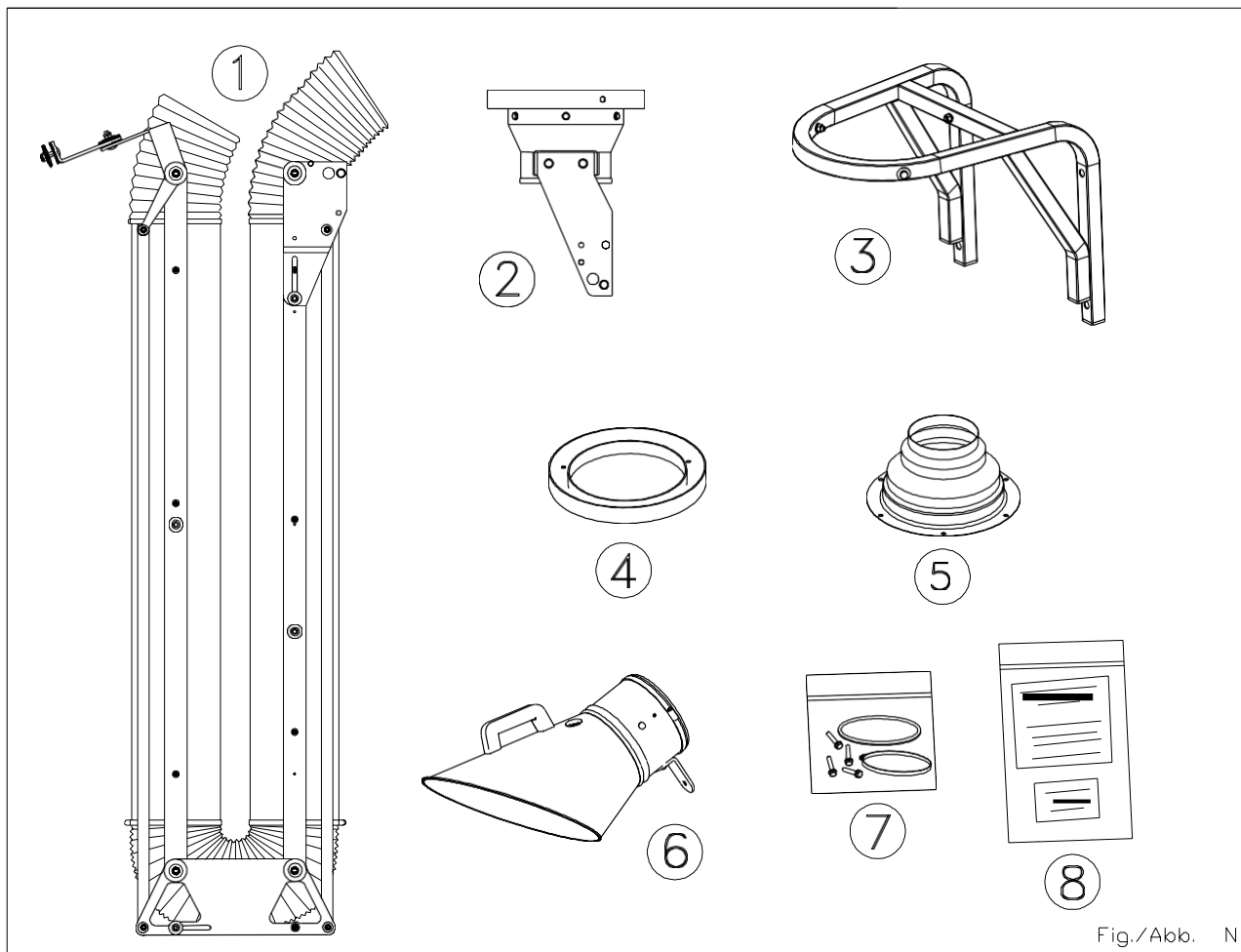
Fig./Abb. N

Fig. N- Suction arms are delivered in three packs:
box n. 1 contains items (1) and (8), box n. 2
contains item (2) and box n. 3 contains the remaining
items (*).

EN

- 1) ARM BODY
- 2) ROTATING BASE
- 3) BRACKET FOR WALL MOUNTING
- 4) COUNTER FLANGE
- 5) FITTING
- 6) METAL HOOD
- 7) BAG WITH INSTALLATION KIT
- 8) BAG ON THE OUTSIDE OF THE PACKAGE WITH MANUAL AND
TICKET WITH ARM ID NUMBER

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PACKAGING
EMBALLAGE
IMBALLAGGIO
VERPACKUNG

PACKAGING

EN

Fig. N shows the components inside the packaging with the weights and dimensions shown in Tab. 1.

An adhesive label indicating the following information is found on the outer part of the packaging:

- Sender, recipient, delivery address, type of goods contained and other information required by the carrier.

-  Before handling the packed product from the delivery vehicle, check that the packaging is entirely intact so as to prevent any components from leaking and/or falling out.
-  When unpacking, move the removed material away from the operating areas so as to eliminate the risk of the operator stumbling over it.
-  Under no circumstances must the end components of the package (cardboard, polyethylene, metal) be disposed of in the environment.
-  It is prescribed for disposal to be implemented by using the regular waste disposal channels according to local regulations.

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